


# Accelerometer

Speed of the Hand

Use GoDirect Accelerometer to determine the maximum speed of a hand moving back and forth.

Setting up:

1. Open Graphical Analysis (or choose New Experiment) and choose Sensor Data Collection. Identify your sensor by its Bluetooth ID and click Connect.
2. Use *only* the Z-axis accelerometer channel. Deselect the Force channel. Under the View Options menu choose Meters. You should see a live readout.
3. Place the sensor at rest on the table on its side, with the z-axis parallel to the surface and the y-axis pointing to the ceiling. Then under the Sensor Actions menu choose Zero. The meter should show a value close to zero.

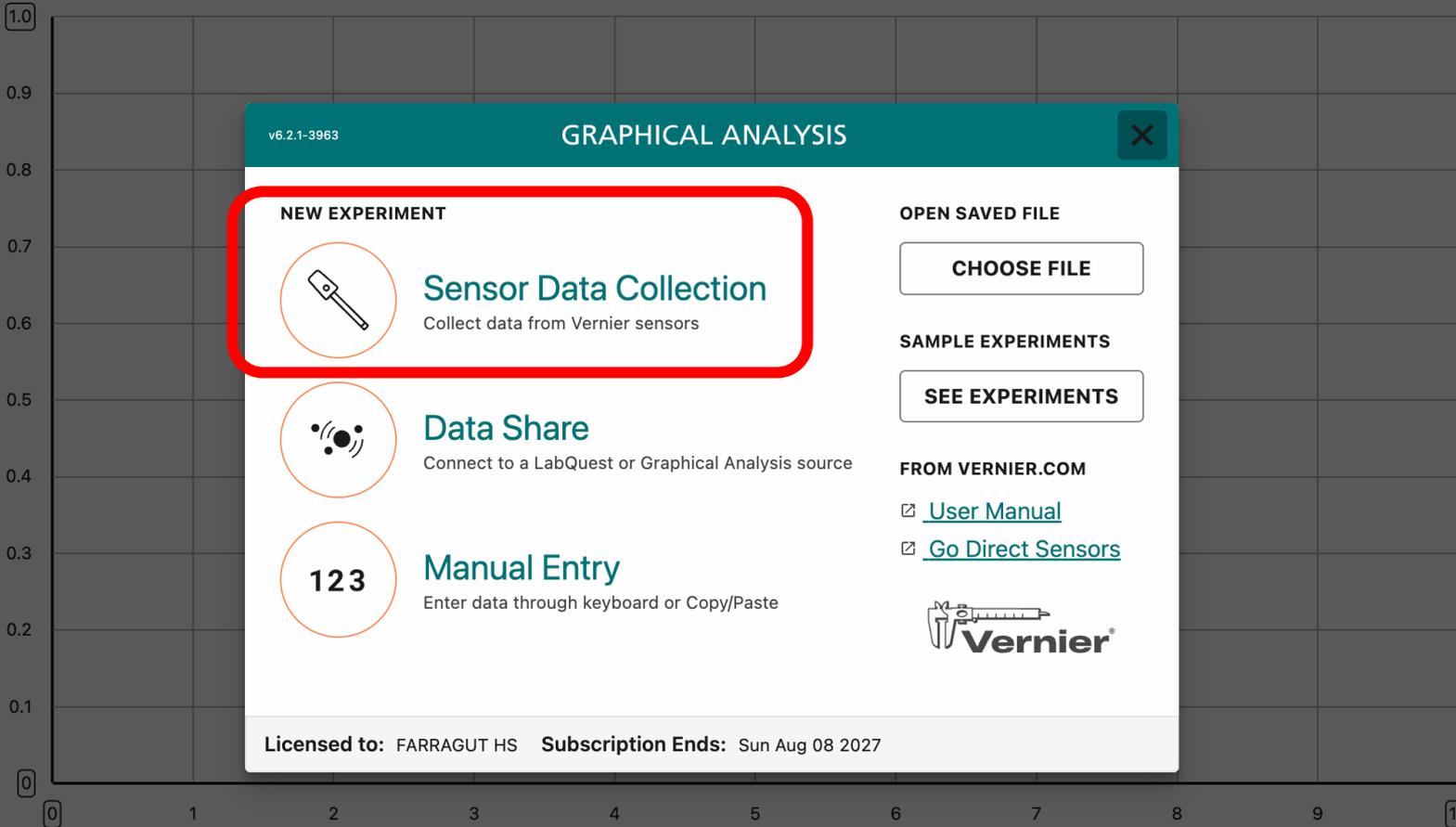

v6.2.1-3963

## GRAPHICAL ANALYSIS

### NEW EXPERIMENT



### Sensor Data Collection

Collect data from Vernier sensors



### Data Share

Connect to a LabQuest or Graphical Analysis source



### Manual Entry

Enter data through keyboard or Copy/Paste

### OPEN SAVED FILE

CHOOSE FILE

### SAMPLE EXPERIMENTS

SEE EXPERIMENTS

### FROM VERNIER.COM


[User Manual](#)

[Go Direct Sensors](#)



Licensed to: FARRAGUT HS    Subscription Ends: Sun Aug 08 2027



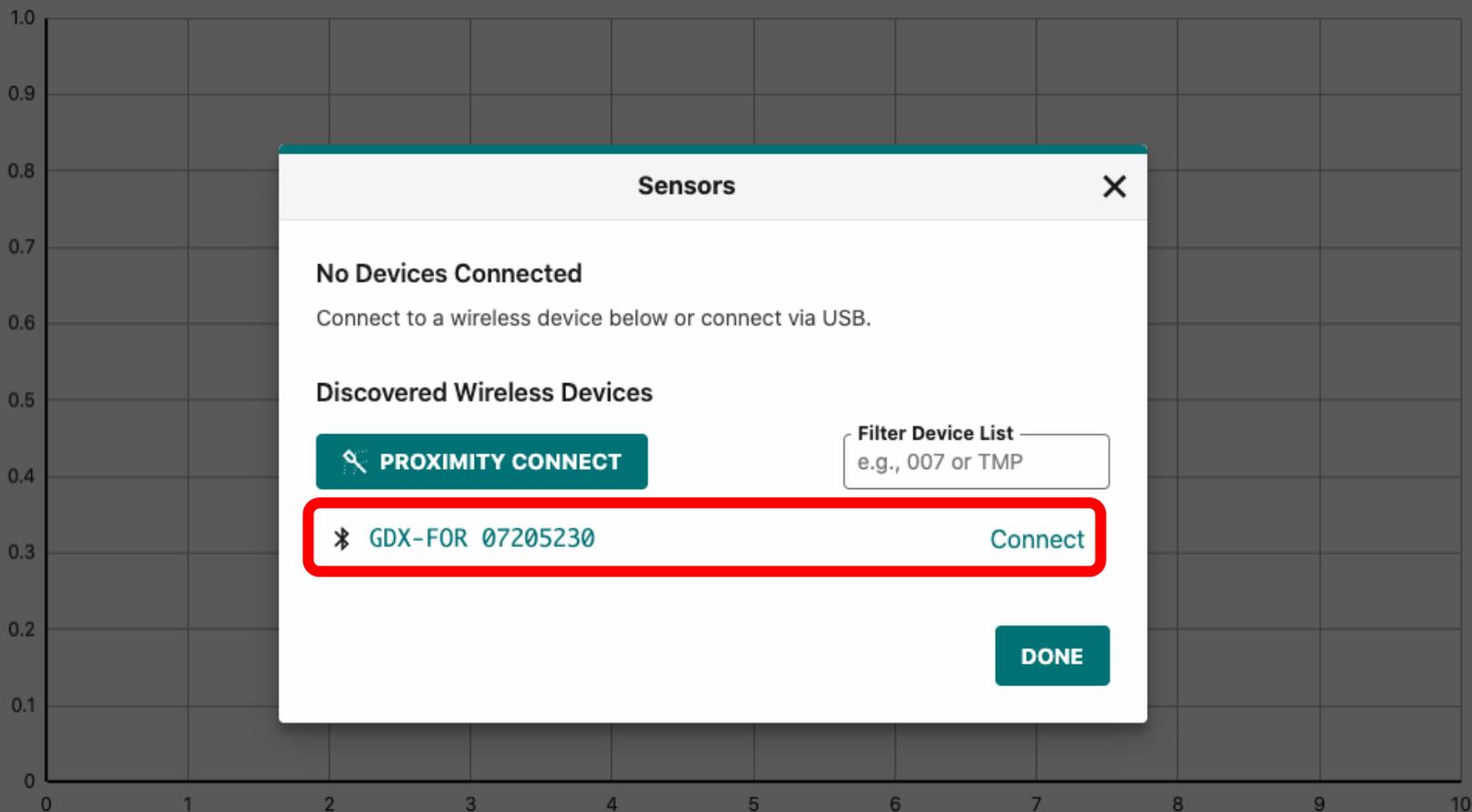

### Sensors



#### No Devices Connected

Connect to a wireless device below or connect via USB.

#### Discovered Wireless Devices

 PROXIMITY CONNECT

Filter Device List

e.g., 007 or TMP

 GDX-FOR 07205230

Connect

DONE





X



Untitled

COLLECT

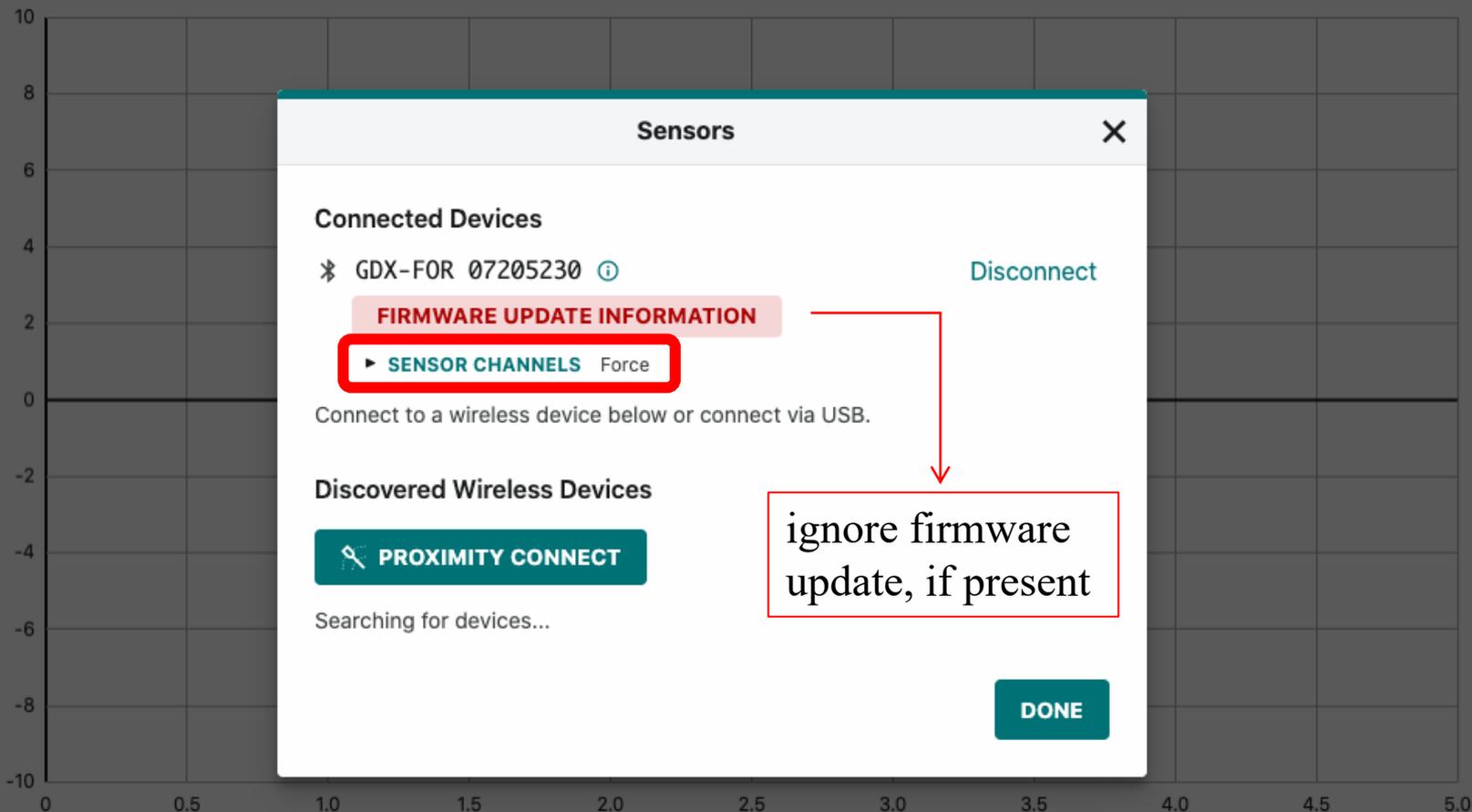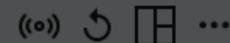

## Sensors



## Connected Devices

\* GDx-FOR 07205230 ⓘ

Disconnect

## FIRMWARE UPDATE INFORMATION

▶ SENSOR CHANNELS Force

Connect to a wireless device below or connect via USB.

## Discovered Wireless Devices

🔍 PROXIMITY CONNECT


Searching for devices...

DONE

ignore firmware  
update, if present


Untitled

COLLECT

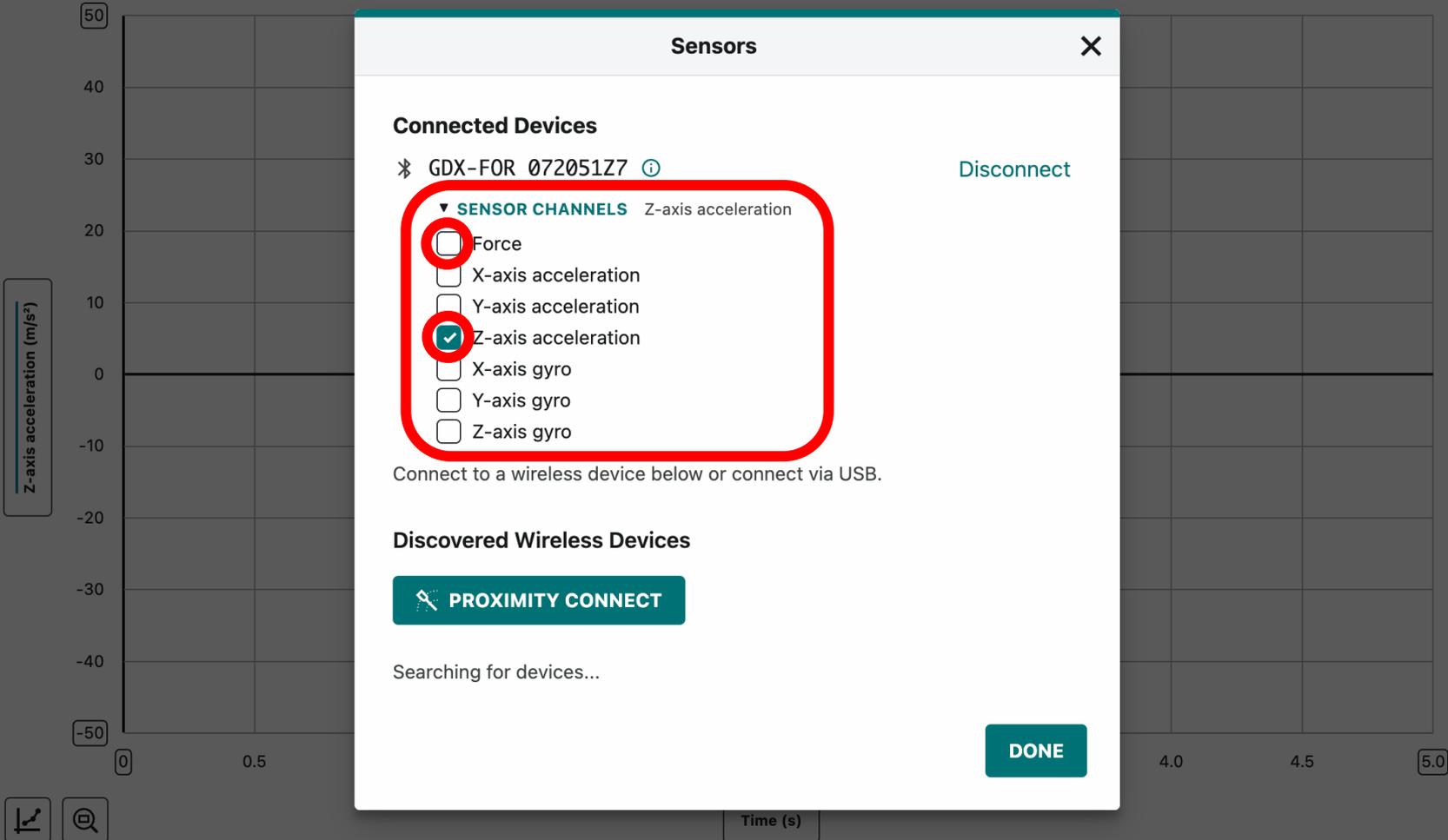



Z-axis acceleration (m/s<sup>2</sup>)

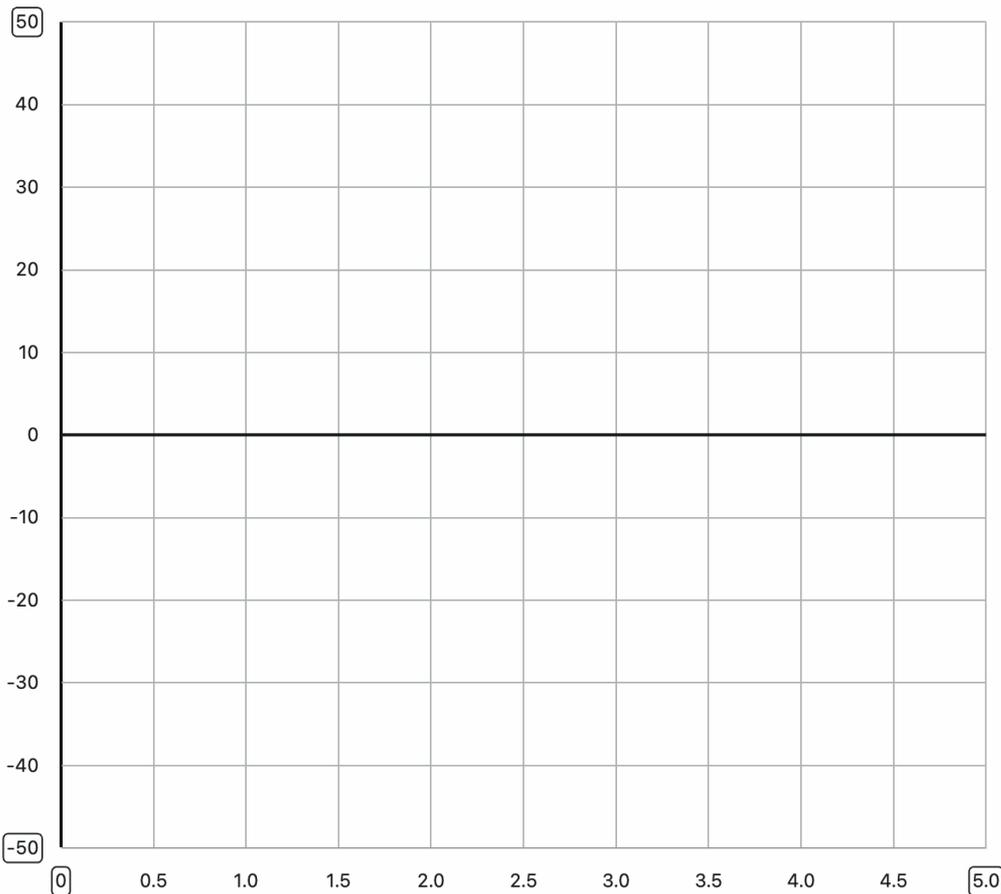

Time (s)

Z-axis accel  
9.68 m/s<sup>2</sup>

- Graph
- 1 Graph ▾
- Data Table
- Meters
- Video
- Notes




Untitled

COLLECT

Help (??) Refresh Home ...

In order to zero the accelerometer, hold the sensor at rest on table as shown here:

Z-axis acceleration: **-0.23 m/s<sup>2</sup>**

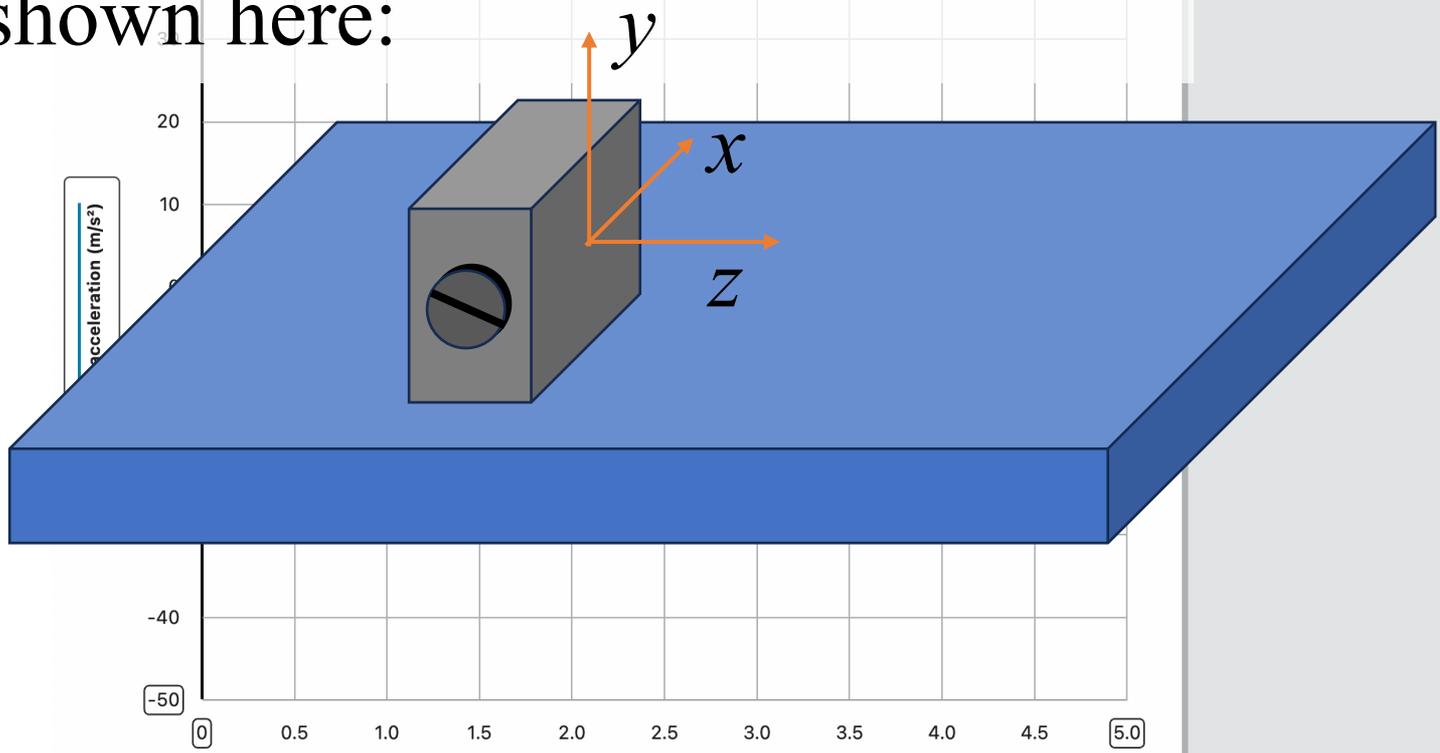

Graph and Zoom icons

Time (s)

Mode: Time Based Rate: 50 samples/s

Z-axis acceleration: **-0.23 m/s<sup>2</sup>**




Untitled

COLLECT


Help (??) Refresh Grid ...

In order to zero the accelerometer, hold the sensor at rest on table as shown here:

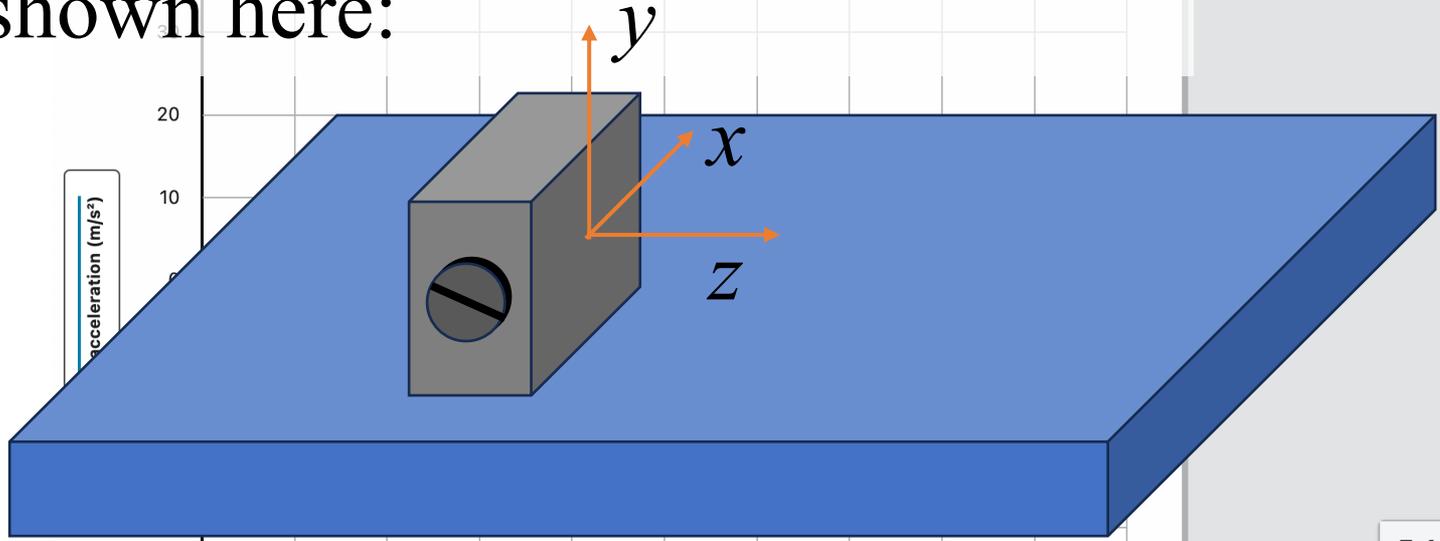

Z-axis acceleration: **-0.15 m/s<sup>2</sup>**

acceleration (m/s<sup>2</sup>)

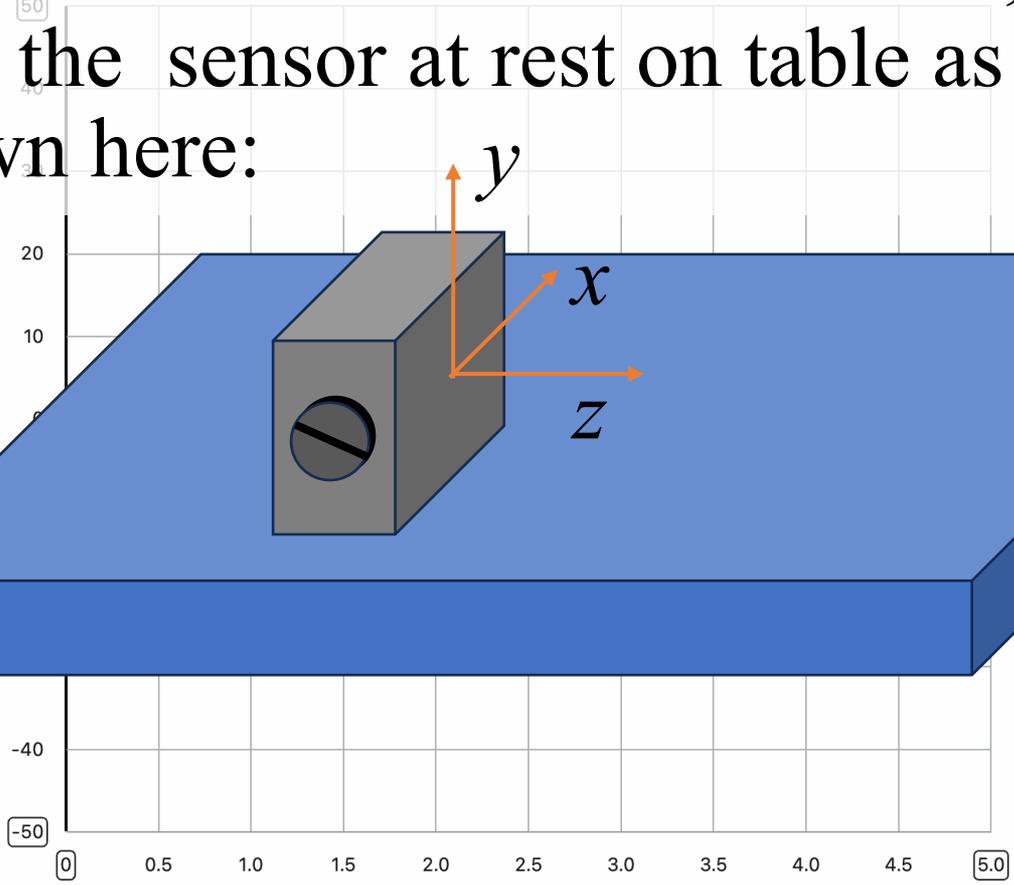

**Z-AXIS ACCELERATION**  
GDX-FOR 072051Z7

Calibrate

**Zero**


Units **m/s<sup>2</sup>**

Z-axis acceleration: **-0.15 m/s<sup>2</sup>**

Graph and Zoom icons

Time (s)

Mode: Time Based Rate: 50 samples/s

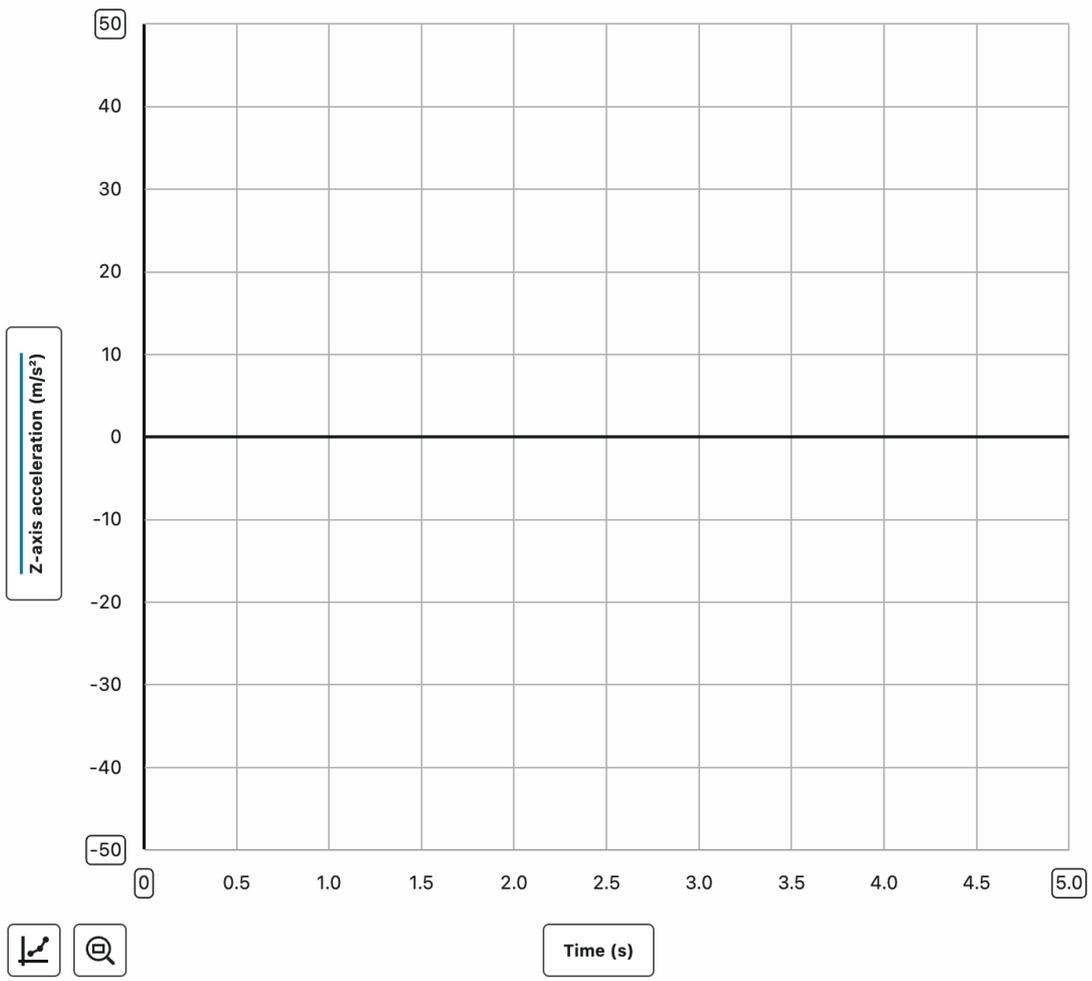

Z-axis acceleration: ...  
**-0.04  $\text{m/s}^2$**



Time (s)





Use GoDirect Accelerometer to determine the maximum speed of a hand moving back and forth.

Doing the experiment:

1. Collect data and move hand quickly back and forth a *distance* of precisely 1.00 m (0.5 m each way). You can move hand back and forth multiple times in one trial.
2. Using the graph and analysis tools determine the *maximum* speed of the hand.
3. Use data from the graph and the *measured distance* to determine the *average* speed of the hand.
4. Based on all available information sketch a graph of velocity vs. time consistent with observations and values.

Click the Collect button and then move the sensor and your hand back and forth several times as shown:

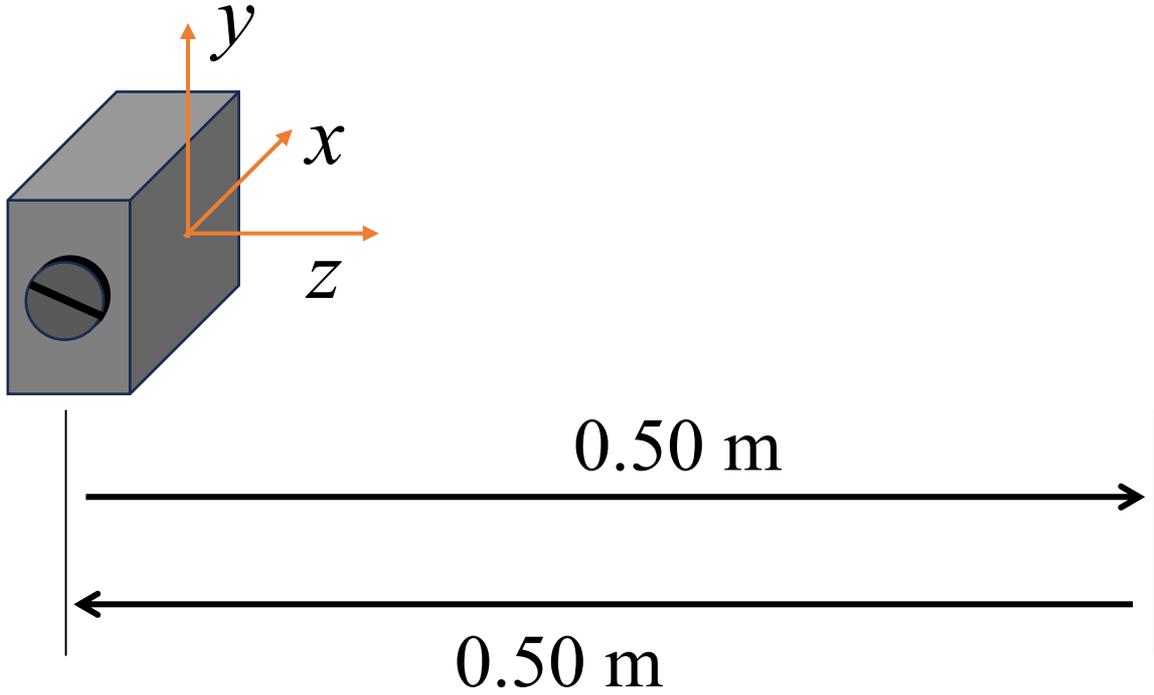

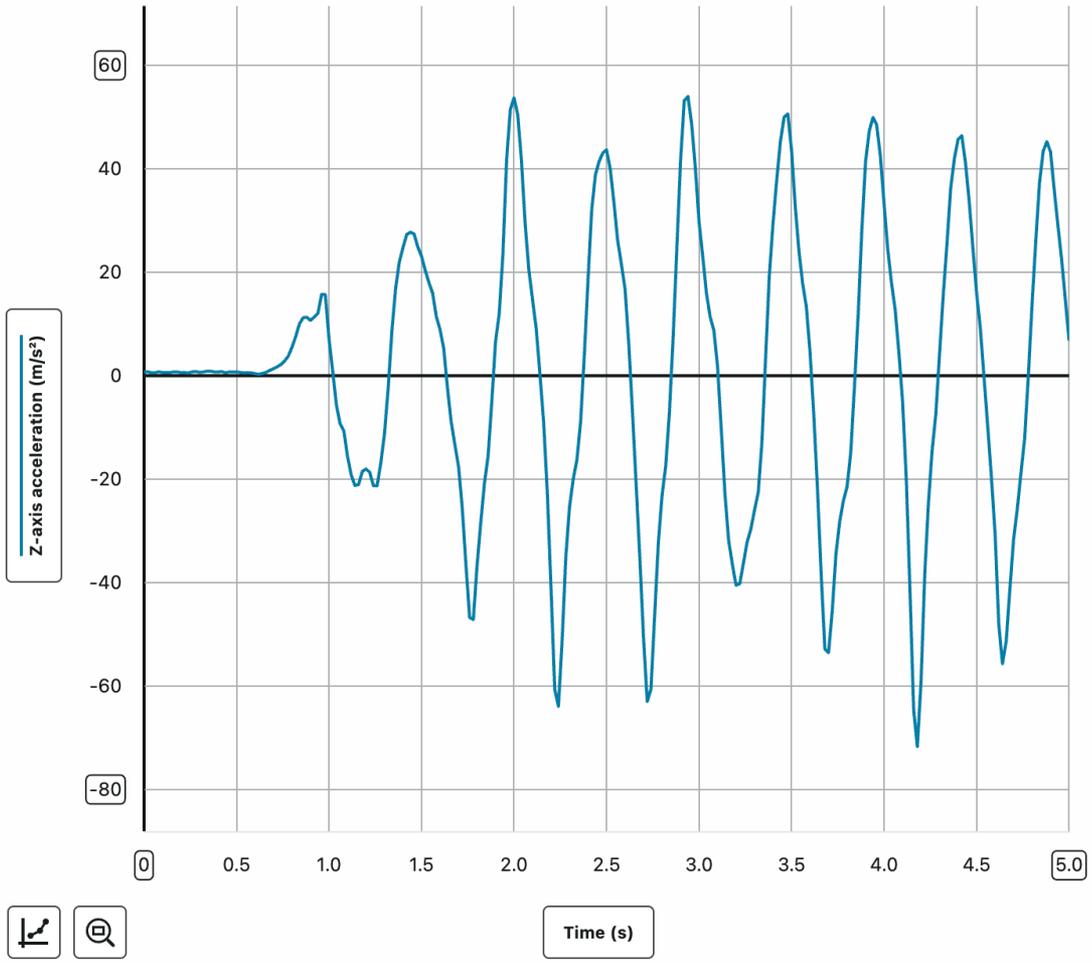

Z-axis acceleration: ...  
**-0.09 m/s<sup>2</sup>**

Typical result should look like this. Repeat as needed until you get some decent, usable data.

Untitled

COLLECT

((o)) ↺ [ ] ...

## Data Collection Settings

✕

Mode Time Based ▾

Time Units s ▾

Rate 50 samples/s

Interval 0.02 s/sample

Start Collection  Manually  
 On a triggering eventEnd Collection  After 5 s duration  
 Manually


Total samples 251

Data Marking ⓘ  Disabled  
 Enabled

CANCEL

DONE

Feel free to  
experiment with  
Data Collection  
Settings

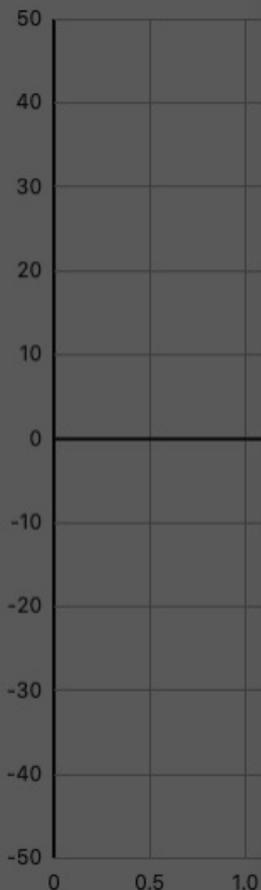Y-axis acceleration (m/s<sup>2</sup>)

Mode: Time Based Rate: 50 samples/s

Y-axis acceleration: -0.02 m/s<sup>2</sup>